

RocketCell™ 3D NSC Xeno-Free Complete Growth Kit



All-in-one ready-to-use kit for 3D neural stem cell culture



Xeno-Free Ready-to-Use, All-in-One Kit

100% synthetic kit with optimal combination of matrix and medium for 2D and 3D NSC culture and maintenance.



Support 3D NSC Network Formation

Promotes 3D neurite and axonal outgrowth for 3D networking formation, ideal for advanced studies.



Easy Cell Harvesting

Simple and efficient cell harvesting with VitroGel® Organoid Recovery Solution

Long-Term Culture and Maintenance




Confidently culture NSCs for over four weeks without affecting their multipotency.

Simple 3D Workflows at Room Temperature



Room temperature operation with simple mixing or coating steps; support lab automation

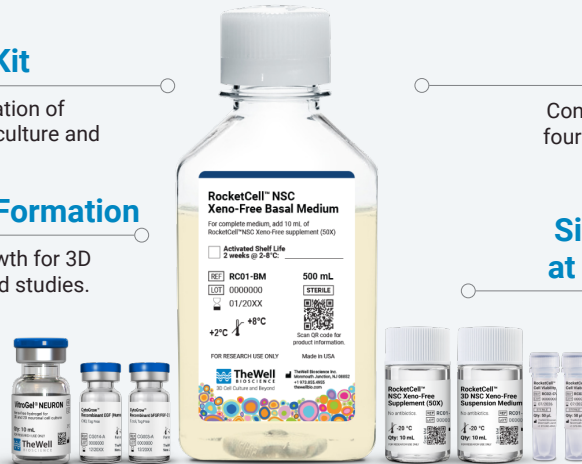

The **RocketCell™ 3D NSC Xeno-Free Complete Growth Kit** is an all-in-one, ready-to-use kit designed to support the expansion and maintenance of neural stem cells (NSCs). This kit offers a fully defined, optimized combination of a synthetic hydrogel and a xeno-free complete medium, eliminating variability and ensuring consistent, reproducible results.

The RocketCell™ 3D NSC Xeno-Free Complete Growth Kit provides a robust, scalable platform for **long-term culture** and **large-scale expansion** of undifferentiated NSCs while maintaining their multipotency and functional integrity. Designed for convenience and performance, this ready-to-use system operates at **room temperature**, integrates seamlessly with **high-resolution imaging** and **high-throughput screening** workflows for advanced neuroscience and regenerative medicine research.

RocketCell™ 3D NSC Xeno-Free Complete Growth Kit (RC01-CGK)

Kit Contents:

VitroGel® NEURON	1 x 10 mL
RocketCell™ 3D NSC Xeno-Free Suspension Medium	1 x 10 mL
RocketCell™ Cell Viability Enhancer (1000X)	2 x 50 µL
RocketCell™ NSC Xeno-Free Basal Medium	1 x 500 mL
RocketCell™ NSC Xeno-Free Supplement (50X)	1 x 10 mL
CytoGrow™ EGF Growth Factor	1 x 10 µg
CytoGrow™ FGF Growth Factor	1 x 10 µg

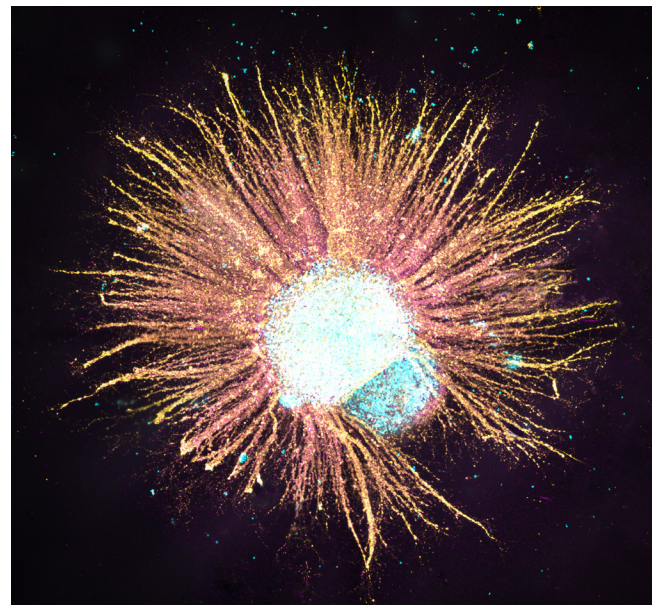


Figure 1. 3D encapsulation culture of NSCs in VitroGel® NEURON hydrogel as NSC spheroids with zonal projections.

Learn more about
**RocketCell™ 3D NSC Xeno-Free
Complete Growth Kit**

thewellbio.com/rocketcell-3d-nsc-xeno-free-complete-growth-kit/



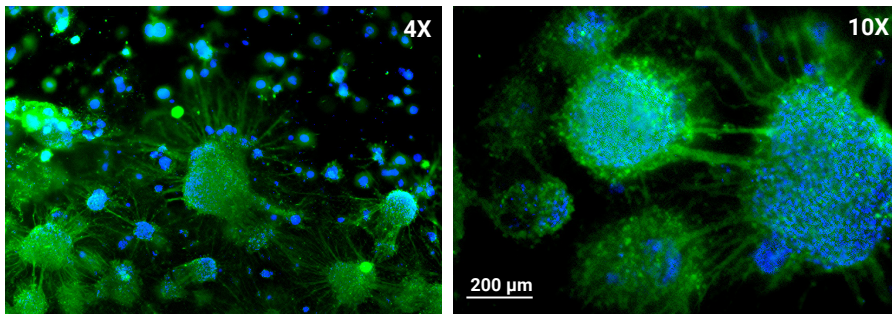


Figure 2. VitroGel® NEURON Hydrogel Supports the 3D Development of Dopaminergic Neurons. iPSCs were cultured in RocketCell™ 3D Xeno-Free IPSC Xeno-Free Growth Kit for 6 passages and were differentiated into Floor Plate Neuro-Epithelial Spheroids.

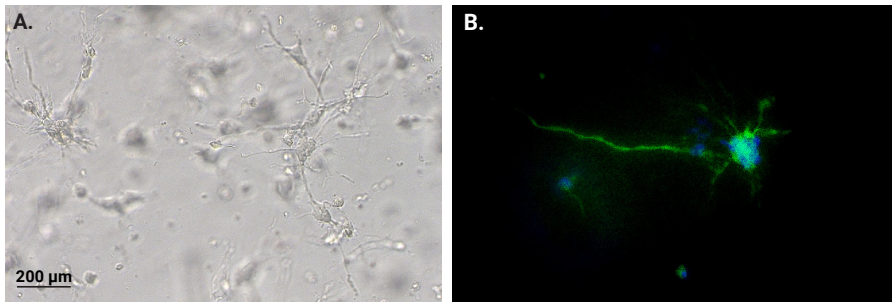


Figure 3. 3D Encapsulation Culture of NSC Cells as Single Cell Suspension in VitroGel® NEURON. VitroGel® NEURON Hydrogel and RocketCell™ NSC Xeno-Free Media support 3D NSC cultures and neurite outgrowth. (A) 3D NSC cultures were established by preparing a single-cell NSC suspension with RocketCell™ NSC Xeno-Free Medium and mixing with VitroGel® NEURON for 6 days. (B) Immunofluorescence staining was performed to evaluate the presence of beta-III-tubulin, a neuron maturation marker. The images were obtained with the Keyence BZ-X microscope at 10 and 20X magnifications.

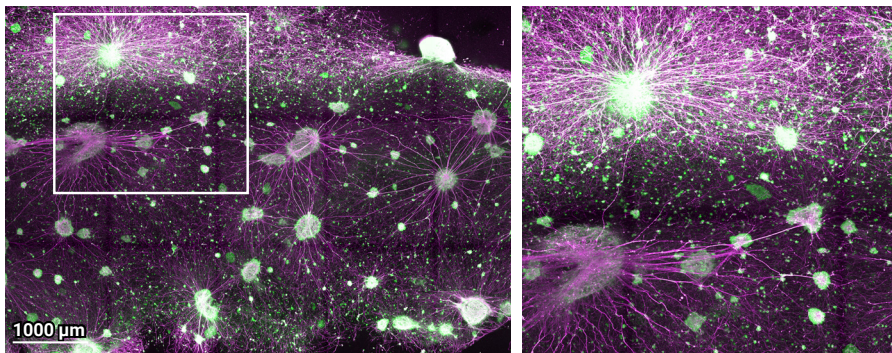


Figure 4. Development of Cholinergic Neurons from iPSCs using FPNE Protocol using VitroGel® NEURON. FPNE Organoids were recovered from VitroGel® STEM using VitroGel® Organoid Recovery Solution and plated with VitroGel® NEURON and RocketCell™ NSC Xeno-Free Growth Medium. Cultures were developed for 14 days prior to indirect immunostaining for Phosphorylated Neurofilament (RT97, purple) and Choline Acetyl Transferase (CHAT, green).

Also Available: **RocketCell™ NSC Xeno-Free Growth Medium (RC01-GM)**

Kit Contents:

RocketCell™ NSC Xeno-Free Basal Medium	1 x 500 mL
RocketCell™ NSC Xeno-Free Supplement (50X)	1 x 10 mL
CytoGrow™ EGF Growth Factor	1 x 10 μg
CytoGrow™ FGF Growth Factor	1 x 10 μg

Related products



RocketCell™ NSC Xeno-Free Growth Medium

Xeno-Free complete medium for 3D and 2D neural stem cell culture



RocketCell™ 3D iPSC Xeno-Free Complete Growth Kit

All-in-one xeno-free kit with optimized matrix, medium, and reagents for 3D iPSC expansion.



RocketCell™ Xeno-Free Supplements

Chemically defined, xeno-free supplements that provide the essential nutrients needed for cell expansion, differentiation, and long-term maintenance.



CytoGrow™ Growth Factors

Premium quality growth factors with research-friendly price for advanced 3D/2D cell culture and beyond.



VitroPrime™ Spread-Attach Plate

Supports exceptional 2D and 3D neuron growth. No PDL-coated plates required.

Plates Available: 6-well, 12-well, 24-well, 48-well, 96-well



VitroPrime™ 3D Culture and Imaging Plate

Premium cover-glass bottom plate for zero disruption 3D cell culture workflow.

Plates Available: 6-well, 24-well, 96-well



VitroGel® Organoid Recovery Solution

Non-enzymatic cell harvesting solution to recover cells or organoids from hydrogel or an animal-based ECM within 15 minutes.



Cyto3D® Live-Dead Assay Kit

Versatile, live/dead cell viability analysis for 3D and 2D cell culture.



Explore our End-to-End Xeno-Free Platform for 3D Cell Culture

VitroGel®
Xeno-Free Hydrogels

CytoGrow™
Premium Growth Factors

RocketCell™
Xeno-Free Culture Media & Supplements

VitroPrime™
Culture Vessels

VitroGel®
Recovery Solution

Cyto3D®
Reagent for 3D Cell Analysis