



3D Cell Culture and Beyond

VitroGel[®] EMT

Low Concentration Hydrogel Protocol for Spheroid, Organoid, and Tumoroid Formation

Synthetic, chemically defined hydrogel that supports spheroid formation, tumoroid culture, and uniquely enables EMT biology

Protocol
Catalog #: VHM08, VHM08S

Revision No: 1.0
Revision Date: 08/2026

www.thewellbio.com
support@thewellbio.com

US TF: +1 866.3D. CELLS
TEL: +1 973.855.4955

For research use only. Not intended for use in diagnostics or therapeutic procedures.



Scan the QR code
to access the latest product updates and resources.

thewellbio.com/product/vitrogel-emt/

TABLE OF CONTENTS

A. Introduction	3
B. Materials	3
C. Gel Preparation	4
D. Xeno-Free Option	4
E. Cell Preparation	5
F. Cell Seeding	5
G. Culture Maintenance	5
H. Downstream Assays	6
Appendix: Related Products	6



VitroGel® EMT

Low-concentration hydrogel protocol for spheroid, organoid, and tumoroid formation



TheWell
BIOSCIENCE

Using VitroPrime™ Ultra-Low Attachment, U-Bottom, 96-Well Plate

A. Introduction

This protocol describes a low-concentration VitroGel® EMT method for forming 3D spheroids, tumoroids, and organoids in the VitroPrime™ Ultra-Low Attachment (ULA), U-Bottom, 96-well plate. By combining a diluted hydrogel scaffold with the gravity-assisted U-bottom geometry, this approach promotes compact cell aggregate formation — even for cell types that do not readily self-aggregate under standard ULA conditions.

Two final gel concentrations are supported:

- 5% VitroGel® EMT — prepared from a 10% intermediate, mixed 1:1 with cells (recommended starting condition)
- 2.5% VitroGel® EMT — prepared from a 5% intermediate, mixed 1:1 with cells (for softer scaffold or faster aggregation)

B. Materials

• TheWell Bioscience Products

- » VitroGel® EMT (Catalog #: VHM08)
- » VitroPrime™ Ultra-Low Attachment, U-Bottom, 96-Well Plate (Catalog #: VP-ULA96U-8)
- » RocketCell™ EMT Supplement (50X), XF (Catalog #: RC05-S50)
(for xeno-free workflows — see Section D)
- » CytoGrow™ Growth Factors (cell-type specific — see Section G)
- » Cyto3D® Live-Dead Assay Kit (Catalog #: BM01) (for downstream viability assessment — see Section H)

• General Lab Supplies

- » Sterile 1.5 mL microcentrifuge tubes
- » Low-retention pipette tips; multichannel or repeater pipette (P200/P300)
- » Centrifuge (300 × g), CO₂ incubator (37°C / 5%), inverted microscope
- » Cell culture medium of choice

C. Gel Preparation

1. Allow VitroGel® EMT to equilibrate to room temperature (5–10 min).
2. Select your target final gel concentration (e.g., 2.5% or 5%) for spheroid formation.
3. Prepare the hydrogel intermediate solution at 2X the final concentration as shown in the table below. The solution will be mixed with cells in a 1:1 ratio (refer to Section F).

Target Final Concentration	Intermediate to Prepare	VitroGel® EMT	Complete Medium
5% (recommended)	10%	25 µL	225 µL
2.5%	5%	12.5 µL	237.5 µL

4. Combine VitroGel® EMT and complete medium in a sterile tube at the ratio above.
5. Triturate 4 times with a small-volume pipette, then switch to a P1000 and triturate 10 more times until homogenous.

NOTE: Prepare the gel intermediate fresh before the experiment. Mix with cells immediately before seeding — do not pre-mix in bulk. Replace the FBS-containing medium with your base medium supplemented with RocketCell™ EMT Supplement (50X), XF (Catalog #: RC05-S50) at the recommended 1X final working concentration.

D. Xeno-Free Option

For fully animal-component-free workflows, replace serum-containing medium with a xeno-free alternative throughout this protocol

XENO-FREE OPTION: Replace FBS-containing medium with your base medium supplemented with RocketCell™ EMT Supplement (50X), XF (Catalog #: RC05-S50) at the recommended working concentration.

Use this xeno-free medium in all gel dilution, cell resuspension, and feeding steps.

Pair with the appropriate CytoGrow™ growth factors for your cell type to support proliferation and phenotype maintenance without serum (see Section G).

RocketCell™ Organoid , AOF Essential-Core Medium is also a compatible base medium for organoid applications.

E. Cell Preparation

1. Harvest cells using your standard dissociation method and quench the enzyme in complete medium.
2. Centrifuge at 300 × g for 3 min, aspirate supernatant.
3. Count cells with a cell viability dye such as Trypan Blue.
4. Resuspend at 2X the intended final seeding density in complete medium.

OPTIONAL: For sensitive cells (e.g., primary cells), add the RocketCell™ Cell Viability Enhancer (1000X) to the medium to achieve a 1X final concentration

F. Cell Seeding

Final seeding volume: 100 μ L per well. Prepare 10-20% extra volume to account for dead volume.

1. Gently mix gel intermediate with cell suspension at 1:1 ratio and pipette up and down 8–10 times. Do not vortex. See suggested volumes in the table below.

Component	Per Well	For 96-well plate (+ 10% extra)
2X Cell Stock	50 μ L	5.28 mL
Gel Intermediate (10% $\rightarrow$ 5% final, or 5% $\rightarrow$ 2.5% final)	50 μ L	5.28 mL
Final volume per well	100 μL	9.6 mL

2. Dispense 100 μ L of the mixture per well of the VitroPrime™ ULA, U-Bottom, 96-well Plate using a multichannel or repeater pipette.
3. Optional: Seal the plate with parafilm and centrifuge at 300 $\times$ g for 3 min to concentrate cells at the well center.
4. Transfer to the incubator at 37°C / 5% CO₂. Do not disturb the cultures for at least 24 hours.

G. Culture Maintenance

Day 1: Inspect under an inverted microscope. Compact aggregation should be visible within 24 hours for most cell types.

- Medium changes: replace 50% of the medium every 2–3 days by removing it from the well periphery with a micropipette. Dispense fresh warm medium along the well wall.
- For defined or xeno-free culture, supplement medium with CytoGrow™ growth factors appropriate for your cell type (e.g., EGF, FGF-2, HGF, VEGF). CytoGrow™ covers 100+ recombinant proteins across all major growth factor families, validated in cancer, stem cell, and organoid models.

H. Downstream Assays

The following TheWell Bioscience products are optimized for in-well analysis of spheroids formed by this protocol:

Product	Application
Cyto3D® Live-Dead Assay Kit	One-step AO/PI viability staining; optimized for 3D spheroids, tumoroids, and organoids; no pre-mixing; add directly to well and image after 5–15 min at 37°C
VitroGel® Cell Recovery Solution	Non-enzymatic dissociation of spheroids from gel matrix for harvest, replating, or downstream analysis

Appendix: Related Products

Product	Catalog #	Description
VitroGel® GBM EMT Kit, XF	VHM08-K	Complete kit for GBM tumoroid formation with xeno-free method
RocketCell™ EMT Supplement (50X), XF	RC05-S50	Xeno-free serum replacement for spheroid and tumoroid formation
RocketCell™ S1 Supplement (50X), Minus Vitamin A, AOF	RCS1-NA	Animal origin-free B27-like Supplement without vitamin A
RocketCell™ S2 Supplement (100X), AOF	RCS2	Animal origin-free N2-like supplement
RocketCell™ S3 Supplement (50X), AOF	RCS3	Animal origin-free N@-B27 like premixed supplement
RocketCell™ Organoid Essential-Core Medium, AOF	RC04-OCM	Universal animal origin-free base medium; replaces N2/B27 for organoid workflows
VitroPrime™ Ultra-Low Attachment, U-Bottom, 96-Well Plate	VP-ULA96U	Ultra-low attachment plate with U-bottom for spheroid formation
CytoGrow™ Growth Factors	CG-	Over 100+ types of premium recombinant proteins across all major growth factor families
VitroGel® Organoid Recovery Solution	MS04-100	Non-enzymatic cell recovery from VitroGel® matrices and animal-based ECMs
Cyto3D® Live-Dead Assay Kit	BM01	Fast, in-well viability staining for 3D spheroids and organoids



Revision No: 1.0
Revision Date: 08/2026

For research use only. Not intended for use in diagnostics or therapeutic procedures.

THEWELL BIOSCIENCE INC. DISCLAIMS ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE. TO THE FULLEST EXTENT ALLOWED BY LAW, IN NO EVENT SHALL THEWELL BIOSCIENCE INC BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF, WHETHER OR NOT FORESEEABLE AND WHETHER OR NOT THEWELL BIOSCIENCE INC. IS ADVISED OF THE POSSIBILITY OF SUCH DAMAGES.

Terms and Conditions: thewellbio.com/terms-and-conditions/

TheWell Bioscience Inc.,
1 Deerpark Dr, Ste C Monmouth Junction, NJ 08852

www.thewellbio.com
support@thewellbio.com

US TF: +1 866.3D. CELLS
TEL: +1 973.855.4955

© 2026 TheWell Bioscience, Inc. All rights reserved. For research use only. Not intended for use in diagnostics or therapeutic procedures. VitroGel®, VitroINk®, Cyto3D®, CytoGrow™, RocketCell™, and VitroPrime™ are trademarks of TheWell Bioscience Inc. All other product names, logos, brands, trademarks, and registered trademarks are property of their respective owners.